



QUICK INFO – SET UP

Fridge-tag 3

1 Packaging content



Fridge-tag



External sensor



Velcro closures (2 pcs)



USB cable (3 m)



USB power supply



Cable channel

Note: Place the external sensor in its predetermined location, before activating the device. See chapter 3 for more information.

Inserting the SIM Card?

See user manual in chapter «[Learn More](#)»
(Only necessary if the SIM card has not been pre-installed at the factory.)

2 State of delivery

You receive the Fridge-tag in sleep mode. The display (LCD) is blank.



Note: If required, note the date of activation on the back of the Fridge-tag.

3 Placing the External Sensor

Two hours before activating the device the external sensor must be positioned in its predetermined location. It is important to place the external sensor in the center of the refrigerator for optimal temperature observation and to avoid any incorrect measurements when starting the device.



4 Activation of the device

Note: Once the device has been activated, operation can neither be interrupted nor paused.

To activate the device, press the **SET** and **READ** buttons simultaneously for more than 3 seconds. The following calendar format will appear: **dd.mm.yyyy**.



If you want to activate the device and the display shows “un def”, the device is not configured.



Need help with initial device configuration?

See user manual in chapter «[Learn More](#)»

5 Setting the calendar format

Press **READ** to change the calendar format.



Option 1

Setting the calendar format to: **dd.mm.yyyy**

Option 2

Setting the calendar format to: **mm.dd.yyyy**

After you have made your selection, press **SET** to save the calendar format.

Note: Depending on the version used, after setting the calendar format, the first digit of the date will start flashing to set date and time.

READ The **READ** button is used to adjust the numbers. Each time you press the **READ** button, the digit that is flashing will increase by 1.

SET The **SET** button is used to save the number. After pressing the **SET** button, the next digit will start flashing.

As soon as the last digit of the time setting is confirmed, the activation is complete.

Need to correct or change settings?

See user manual in chapter «[Learn More](#)»

6 Activation completed




The activated Fridge-tag must be placed immediately in its predetermined location. Connect the device with the external sensor. For max. 1 minute after activation no temperature is displayed on the screen.

Connect included USB cable with the Fridge-tag and the power supply.

Placing the Fridge-tag or connection error?

See user manual in chapter «[Learn More](#)»

7 Fridge-tag information

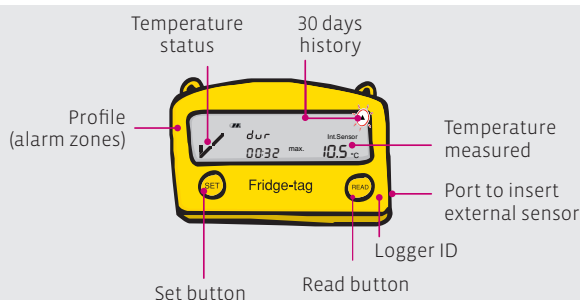

Learn More

Setting date and time



Device configuration



Insert SIM Card



Change settings



Connection error



Placing the Fridge-tag



Scan the QR code or click on www.berlinger.com/support/fridge-tag3

& berlinger


Berlinger & Co. AG
Switzerland
www.berlinger.com



1 Temperature status


Indication of an OK display

In measurement mode, when no alarm has been triggered, the following information is displayed:

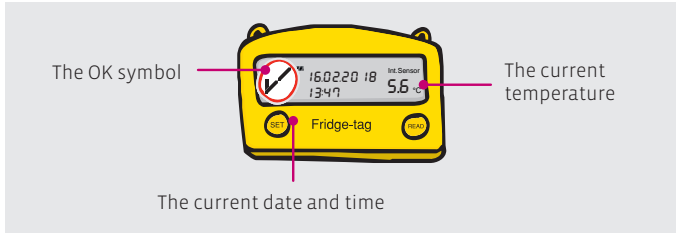


Indication of an alarm display

In measurement mode, when an alarm has been triggered, the following information is displayed:

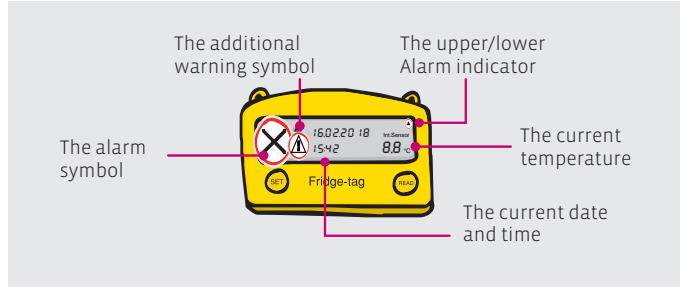


2 Reading the history

Example of an OK display

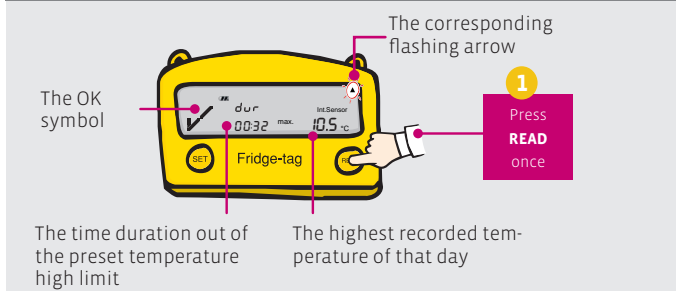


Example of an alarm display

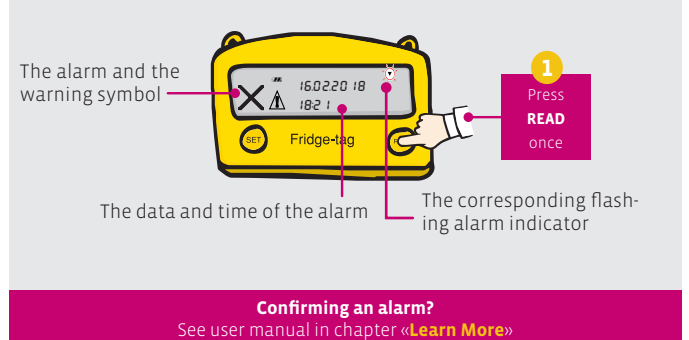

Confirming an alarm?
See user manual in chapter «Learn More»

Alarm display: additional information

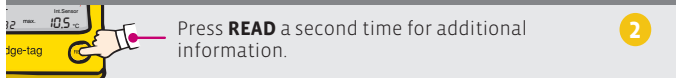


Note: Continue repetitively pressing the **READ** button to read out the details of the past 30 days.



Note: Press and hold the **READ** button again for at least 3 seconds and the next alarm event will appear on the screen.

Need help with Readout mode?
See user manual in chapter «Learn More»

3 Download reports

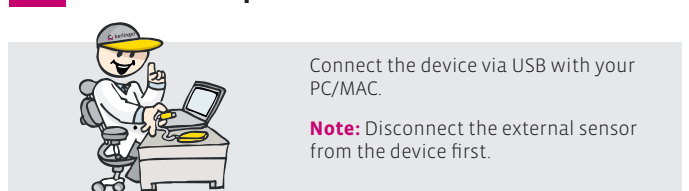


Note: Disconnect the external sensor from the device first.

The Fridge-tag will now generate a PDF and an ASCII report. This process can take up to 2 minutes. After download, please connect the device with the external sensor again.

Looking to download reports?
See user manual in chapter «Learn More»


4 SMS status query

Sending an SMS to the device with the message "status" triggers a response with the status OK or Alarm and the current temperature.

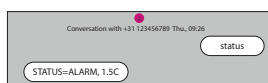

5 Alarm confirmation

Reading out the alarm on the device generates a confirmation notification.

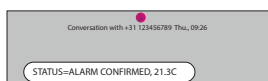

Setting options via SMS?
See user manual in chapter «Learn More»

6 Cloud Version

berlinger smartview Log in to your Data Management Software (Berlinger Smart-View etc.) to check the configuration status and the data.

Learn More

<p>Confirming an alarm</p>	<p>Readout mode</p>	<p>Download report</p>	<p>SMS setting options</p>	<p>Fridge-tag support</p>	<p>Product information</p>
-----------------------------------	----------------------------	-------------------------------	-----------------------------------	----------------------------------	-----------------------------------