HANDLING INSTRUCTION

berlinger® lausanne KIT Universal Opener



Berlinger Special AG

Mitteldorfstrasse 2 9608 Ganterschwil

Tel. +41 71 982 88 11 info.special@berlinger.comwww.berlinger.com



HANDLING INSTRUCTION

WHAT IS INCLUDED

In this Universal Opener Berlinger Lausanne Kit box, you will find:

• One (1) Berlinger Lausanne Kit Universal Opener, one (1) Power cable, one (1) Lausanne test kit, one (1) slip-resistent mat.

Note: total weight approx. 32.0 kg



PREPARATION

Please make sure that:

- The content of the Universal Opener Berlinger Lausanne Kit is complete.
- 1. Connect the machine with the air pressure. Minimum needed pressure is 4 bar.
- 2. Connect the machine to the electricity by plugging the cable. The standard power range is from 100 to 240 VAC (50/60 Hz).
- 3. Use the two handles on each side of the opener to lift it.



HANDLING THE UNIVERSAL OPENER BERLINGER LAUSANNE-KIT

- 4. Turn the power on by pushing the switch on the back and check first, if the container cap is firmly and properly closed.
- 5. Place the Berlinger Lausanne Kit firmly into the white cup.

Note: Do not grab into the machine, sharp cutting wheels inside.

- 6. To start the opening process slide down the door by using the black leveron the front. All is protected by the door and you must take away your hands. After 6-8 seconds the process is finished and the door opens automatically.
- 7. The door is open and the Berlinger Lausanne-Kit ready to be taken out.
- 8. Open the cap by turning it in an anti-clockwise direction.
- 9. The separated top of the cap can be re-used to close the container again.

Note: Protect your hands by opening the cap with the slip-resistant mat.











WARNING

- Keep your hands away from places that bear this warning sign.
 There is a risk that hands may be injured.
- · Use eye protection.



www.berlinger.com/handling-instructions

For additional information, check out the Berlinger Lausanne Kit product webpage

www.berlinger.com/berlinger-lausanne-kit

Contact us for the detailed handling instruction or user manual to replace the cutting wheels.

info.special@berlinger.com

